

SuperScript™ II Reverse Transcriptase

Cat. No. 18064-022

Size: 2,000 units

Cat. No. 18064-014

Size: 10,000 units

Cat. No. 18064-071

Size: 4 × 10,000 units

Conc. 200 U/μL

Store at -20°C (non-frost-free)

Description

SuperScript™ II Reverse Transcriptase (RT) is an engineered version of M-MLV RT with reduced RNase H activity and increased thermal stability. The enzyme is purified to near homogeneity from *E. coli* containing the modified *pol* gene of Moloney Murine Leukemia Virus (1,2). The enzyme can be used to synthesize first-strand cDNA at higher temperatures than conventional M-MLV RT, providing increased specificity, higher yields of cDNA, and more full-length product. It can generate cDNA up to 12.3 kb.

Components

SuperScript™ II RT, 5X First-Strand Buffer (250 mM Tris-HCl, pH 8.3 at room temperature; 375 mM KCl; 15 mM MgCl₂), 0.1 M DTT

Storage Buffer

20 mM Tris-HCl (pH 7.5), 100 mM NaCl, 0.1 mM EDTA, 1 mM DTT, 0.01% (v/v) NP-40, 50% (v/v) glycerol

Storage Conditions

Store all components at -20°C in a non-frost-free freezer.

Thaw 5X First-Strand Buffer and 0.1 M DTT at room temperature just prior to use and refreeze immediately.

Unit Definition

One unit incorporates 1 nmole of dTTP into acid-precipitable material in 10 min. at 37°C using poly(A)•oligo(dT)₂₅ as template-primer (3).

Intended Use

For research use only. Not intended for any animal or human therapeutic or diagnostic use.

Part no. 18064.pps

MAN0001342

Rev. Date: 20 May 2010

First-Strand cDNA Synthesis Using SuperScript™ II RT

A 20- μ L reaction volume can be used for 1 ng–5 μ g of total RNA or 1–500 ng of mRNA.

- Add the following components to a nuclease-free microcentrifuge tube:

Oligo(dT) ₁₂₋₁₈ (500 μ g/mL) <i>or</i>	1 μ L
50–250 ng random primers <i>or</i>	
2 pmole gene-specific primer (GSP)	
1 ng to 5 μ g total RNA <i>or</i>	x μ L
1–500 ng of mRNA	
1 μ L dNTP Mix (10 mM each)	1 μ L
Sterile, distilled water	to 12 μ L
- Heat mixture to 65°C for 5 min and quick chill on ice. Collect the contents of the tube by brief centrifugation and add:

5X First-Strand Buffer	4 μ L
0.1 M DTT	2 μ L
RNaseOUT™ (40 units/ μ L) (optional)*	1 μ L

*RNaseOUT™ (Cat. No. 10777-019) is required if using <50 ng starting RNA.
- Mix contents of the tube gently. If you are using oligo(dT)₁₂₋₁₈ or GSP, incubate at 42°C for 2 min. If you are using random primers, incubate at 25°C for 2 min.
- Add 1 μ L (200 units) of SuperScript™ II RT and mix by pipetting gently up and down.
If you are using less than 1 ng of RNA, reduce the amount of SuperScript™ II RT to 0.25 μ L (50 units) and add sterile, distilled water to a 20 μ L final volume.
If you are using random primers, incubate tube at 25°C for 10 min.
- Incubate at 42°C for 50 min.
- Inactivate the reaction by heating at 70°C for 15 min.

First-Strand cDNA Synthesis Using SuperScript™ II RT, Continued

The cDNA can now be used as a template for amplification in PCR. However, amplification of some PCR targets (>1 kb) may require the removal of RNA complementary to the cDNA. To remove RNA complementary to the cDNA, add 1 μL (2 units) of *E. coli* RNase H and incubate at 37°C for 20 min.

PCR

The following is intended as a guideline and starting point when using first-strand cDNA in PCR with *Taq* DNA polymerase. The optimal concentration of Mg^{++} will vary depending on the template and primer pair.

Use only 10% of the first-strand reaction for PCR. Higher volumes may not increase amplification and may result in decreased amounts of PCR product.

1. Add the following to a PCR tube:

10X PCR Buffer [200 mM Tris-HCl (pH 8.4), 500 mM KCl]	5 μL
50 mM MgCl_2	1.5 μL
10 mM dNTP Mix	1 μL
Forward primer (10 μM)	1 μL
Reverse primer (10 μM)	1 μL
<i>Taq</i> DNA polymerase (5 U/ μL)	0.4 μL
cDNA from first-strand reaction	2 μL
autoclaved, distilled water	to 50 μL

- Mix gently and layer with 1–2 drops (~50 μL) of silicone oil. (*Note: silicone oil is unnecessary in thermal cyclers equipped with a heated lid.*)
- Heat reaction to 94°C for 2 min to denature.
- Perform 15 to 40 cycles of PCR. Use the recommended annealing and extension conditions for your *Taq* DNA polymerase.

Product Qualification

The Certificate of Analysis provides detailed quality control information for each product. Certificates of Analysis are available at www.invitrogen.com/support.

Additional Products

RNaseOUT™ Recombinant Ribonuclease Inhibitor (40 units/ μ L) is available separately from Invitrogen (Cat. no. 10777-019).

References

1. Kotewicz, M.L., D'Alessio, J.M., Driftmier, K.M., Blodgett, K.P., and Gerard, G.F. (1985) *Gene* 35, 249.
2. Gerard, G.F., D'Alessio, J.M., Kotewicz, M.L., and Noon, M.C. (1986) *DNA* 5, 271.
3. Houts, G.E., Miyagi, M., Ellis, C., Beard, A., and Beard, J.W. (1979) *J. Virol.* 29, 517.
4. Kotewicz, M.L., Sampson, C.M., D'Alessio, J.M., and Gerard, G.F. (1988) *Nuc. Acids Res.* 16, 265.

Limited Use Label Licenses

Limited Use Label License No. 5: Invitrogen Technology

The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) not to transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. For products that are subject to multiple limited use label licenses, the terms of the most restrictive limited use label license shall control. Life Technologies Corporation will not assert a claim against the buyer of infringement of patents owned or controlled by Life Technologies Corporation which cover this product based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Life Technologies is willing to accept return of the product with a full refund. For information about purchasing a license to use this product or the technology embedded in it for any use other than for research use please contact Out Licensing, Life Technologies, 5791 Van Allen Way, Carlsbad, California 92008 ; Phone (760) 603-7200 or e-mail: outlicensing@lifetech.com

©2010 Life Technologies Corporation. All rights reserved.