

## M-MLV Reverse Transcriptase

**Cat. Nos.**

28025-013

28025-021

**Size:**

40,000 units

200,000 units

**Conc: 200 U/μl****Store at -20°C (non-frost-free)**

### Description

Moloney Murine Leukemia Virus Reverse Transcriptase (M-MLV RT) uses single-stranded RNA or DNA in the presence of a primer to synthesize a complementary DNA strand. This enzyme is isolated (1) from *E. coli* expressing a portion of the *pol* gene of M-MLV on a plasmid (2, 3). The enzyme is used to synthesize first-strand cDNA up to 7 kb.

### Components

M-MLV RT

5X First-Strand Buffer [250 mM Tris-HCl (pH 8.3 at room temperature),  
375 mM KCl, 15 mM MgCl<sub>2</sub>]

0.1 M DTT

### Storage Buffer

20 mM Tris-HCl (pH 7.5), 100 mM NaCl, 0.1 mM EDTA, 1 mM DTT,  
0.01% (v/v) NP-40, 50% (v/v) glycerol

### Storage Conditions

Store all components at -20°C in a non-frost-free freezer.

### Additional Components

Thaw 5X First-Strand Buffer and 0.1 M DTT at room temperature just prior to use and refreeze immediately.

RNaseOUT™ Recombinant Ribonuclease Inhibitor (40 units/μl) is available separately from Invitrogen (Cat. No.: 10777-019).

### First-Strand cDNA Synthesis Using M-MLV RT

A 20- $\mu$ l reaction volume can be used for 1 ng–5  $\mu$ g of total RNA or 1–500 ng of mRNA.

1. Add the following components to a nuclease-free microcentrifuge tube:
  - 1  $\mu$ l oligo (dT)<sub>12-18</sub> (500  $\mu$ g/ml), or 50–250 ng random primers, or  
2 pmole gene-specific primer
  - 1 ng to 5  $\mu$ g total RNA or 1 ng to 500 ng of mRNA
  - 1  $\mu$ l 10 mM dNTP Mix (10 mM each dATP, dGTP, dCTP and dTTP at neutral pH)
  - Sterile, distilled water to 12  $\mu$ l
2. Heat mixture to 65°C for 5 min and quick chill on ice. Collect the contents of the tube by brief centrifugation and add:
  - 4  $\mu$ l 5X First-Strand Buffer
  - 2  $\mu$ l 0.1 M DTT
  - 1  $\mu$ l RNaseOUT™ Recombinant Ribonuclease Inhibitor (40 units/ $\mu$ l)  
(**Note:** When using less than 50 ng of starting RNA, the addition of RNaseOUT™ is essential.)
3. Mix contents of the tube gently and incubate at 37°C for 2 min.
4. Add 1  $\mu$ l (200 units) of M-MLV RT,<sup>a</sup> and mix by pipetting gently up and down. If using random primers, incubate tube at 25°C for 10 min.
5. Incubate 50 min at 37°C.
6. Inactivate the reaction by heating at 70°C for 15 min.

The cDNA can now be used as a template for amplification in PCR. However, amplification of some PCR targets (>1 kb) may require the removal of RNA complementary to the cDNA. To remove RNA complementary to the cDNA, add 1  $\mu$ l (2 units) of *E. coli* RNase H and incubate at 37°C for 20 min.

<sup>a</sup>If less than 1 ng of RNA is used, reduce the amount of M-MLV RT in the reaction to 0.25  $\mu$ l (50 units), and add the sterile, distilled water to 20- $\mu$ l final volume.

## PCR Reaction

Use only 10% of the first-strand reaction (2  $\mu$ l of the reaction from the previous page) for PCR. Adding larger amounts of the first-strand reaction may not increase amplification and may result in decreased amounts of PCR product.

1. Add the following to a PCR reaction tube for a final reaction volume of 50  $\mu$ l:

- 5  $\mu$ l 10X PCR Buffer [200 mM Tris-HCl (pH 8.4), 500 mM KCl]
- 1.5  $\mu$ l 50 mM  $MgCl_2$ <sup>b</sup>
- 1  $\mu$ l 10 mM dNTP Mix
- 1  $\mu$ l amplification primer 1 (10  $\mu$ M)
- 1  $\mu$ l amplification primer 2 (10  $\mu$ M)
- 0.4  $\mu$ l *Taq* DNA polymerase (5 U/ $\mu$ l)
- 2  $\mu$ l cDNA (from first-strand reaction)
- 38.1  $\mu$ l autoclaved, distilled water

2. Mix gently and layer 1–2 drops (~50  $\mu$ l) of silicone oil over the reaction. (*Note: the addition of silicone oil is unnecessary in thermal cyclers equipped with a heated lid.*)
3. Heat reaction to 94°C for 2 min to denature.
4. Perform 15 to 40 cycles of PCR. Annealing and extension conditions are primer and template dependent and must be determined empirically.

<sup>b</sup> For best results, determine the optimal concentration of  $MgCl_2$  empirically for each template-primer pair.

## Quality Control

This product has passed the following quality control assays: SDS-polyacrylamide gel analysis for purity; functional absence of endodeoxyribonuclease, 3' and 5' exodeoxyribonuclease, and ribonuclease activities; yield and length of cDNA product.

### Unit Definition

One unit incorporates 1 nmole of dTTP into acid-precipitable material in 10 min at 37°C using poly(A)•oligo(dT)<sub>25</sub> as template-primer (4).

### References

1. D'Alessio, J. M., and Gerard, G. F. (unpublished).
2. Kotewicz, M., D'Alessio, J., Driftmeier, K., Blodgett, K., and Gerard, G. (1985) *Gene* 35, 249.
3. Gerard, G. F., D'Alessio, J. M., Kotewicz, M. L., and Noon, M. C. (1986) *DNA* 5:4, 271.
4. Houts, G. E., Miyagi, M., Ellis, C., Beard, A., and Beard, J. W. (1979) *J. Virol.* 29, 517.

### Limited Use Label License No. 4: Products for PCR which do not include any rights to perform PCR

This product is optimized for use in the Polymerase Chain Reaction (PCR) covered by patents owned by Roche Molecular Systems, Inc. and F. Hoffmann-La Roche, Ltd. ("Roche"). No license under these patents to use the PCR process is conveyed expressly or by implication to the purchaser by the purchase of this product. A license to use the PCR process for certain research and development activities accompanies the purchase of certain reagents from licensed suppliers such as Invitrogen, when used in conjunction with an Authorized Thermal Cycler, or is available from Applied Biosystems. Further information on purchasing licenses to practice the PCR process may be obtained by contacting the Director of Licensing at Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404 or at Roche Molecular Systems, Inc., 1145 Atlantic Avenue, Alameda, California 94501.